



LabGadget

www.labgadget.com

x-tracta

The disposable gel extraction tool


User Instructions

This x-tracta tool by LabGadget provides a faster, safer and more convenient way to extract and dispense gel samples.


Attention: Do not squeeze the bulb until dispensing at step 4.

To Dispense: Best to use a quick snap-action squeeze.


1. Insert tool into gel

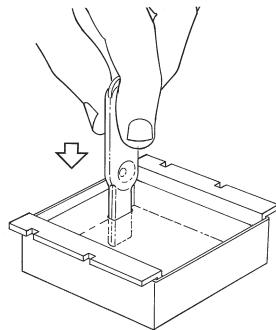


2. Rock slightly to cut into gel

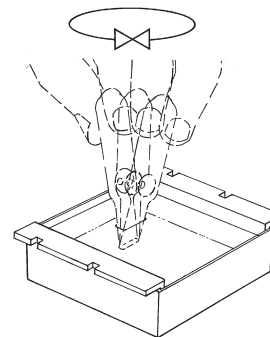

3. Remove Tool at a slant

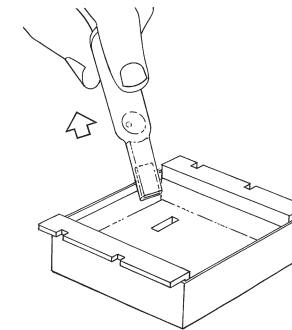

4. Dispense

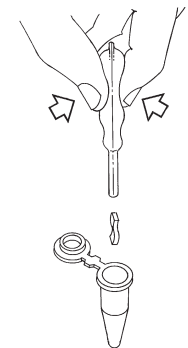

In accordance with good laboratory practice, the x-tracta tool should be used once and discarded to prevent cross-contamination.

LabGadget warrants that its x-tracta tool is free from defects at the time of shipment. LabGadget shall not be liable for any loss of use or profits or facilities or any other indirect, incidental or consequential damages arising out of the use or performance of the product involved.

Patent pending